



Afla- γ [®] AQUA™
QUANTITATIVE STRIP TESTS

INSTRUCTION GUIDE FOR PEANUTS



VICAM[®]

A Waters Business



CULTIVATING SUCCESS THROUGH SCIENCE™

For more than 20 years, VICAM has been the global provider of choice for next-generation food safety technology and rapid mycotoxin test solutions.

The goal of securing the world's food supply requires knowledge, cooperation, and collaboration. To advance that objective, VICAM works closely with government regulatory agencies, international standards bodies, major food industry laboratories, and leading research institutions around the world. With a global scientific and distribution presence, VICAM has access to the latest research advances, industry expertise, and regulatory information needed to design solutions that maximize food safety and quality at every stage of the global food supply chain.

As the trusted partner of the agricultural industry, VICAM is proud to consistently deliver the superior products, responsive service, and individualized support for your industry-specific goals.



AFLA-V AQUA PEANUT KIT 176003776

Kit Contents

Afla-V Strip Tests	100000247
Micro-Pipette Tips, 100 μ L	600001109
Filter Paper	600001522
Extraction Tubes, 70 mL	600001545
Afla-V AQUA Peanut and Peanut Paste Barcodes	715005111

REQUIRED BUT NOT INCLUDED

AQUA Premix Solution (900 mL).....	100000339
AQUA Premix Solution (4 L)	100000345
AQUA Premix Solution (20 L).....	100000347
OR	
AQUA Solution A	100000336
AQUA Solution B	100000337

Vertu Mycotoxin BEQ, 110V

176002078

Vertu Mycotoxin BEQ, 220V

176002079

EQUIPMENT PACKAGE CONTENTS

Vertu Reader, 110V & 220V	725000574
Vertu Bar-Code Scanner	725000576
Printer, 110V & 220V	725000577
Vortex Mixer	23040
Digital Timer	G4036
Micro-Pipettor, 100 μ L.....	600001108
Filter Funnel, 65 mm (4 pack)	36020
Strip Test Vial Rack	600001107
Graduated Cylinder, 50 mL.....	20050
Vertu Basic Instruction Guide.....	715002426

STORAGE AND SAMPLING



STORAGE RECOMMENDATIONS

Strip tests must be refrigerated. Store at 4 ° – 8 °C (39 ° – 46 °F).

Strip test and sample extract should be at room temperature of between 20 °C (68 °F) and 25 °C (77 °F) before use. After removing strips from refrigerator, leave strips at room temperature for 20 minutes or longer to equilibrate.

Use in accordance with VICAM's testing procedure. To be used for testing food products only. Not to be ingested or applied to the skin.

SAMPLING

Mycotoxins tend to occur in scattered locations in large peanut and tree nut loads and may be detectable in only a small percentage of nuts in a lot. The uneven distribution of contaminated nuts can cause test results to vary significantly from sample to sample. It is therefore essential to take a representative sample from the lot. Product must be collected from different locations in a static lot based on a probing pattern. The probe must draw from the top to the bottom of the lot. The samples obtained from the probes must be ground and mixed well and a subsample taken for testing. Contact your local regulatory authorities or VICAM for more information on sampling.



The information contained in each barcode includes the test name, test lot number, method of scanning, parameters of measure, and algorithm for result calculation. This information is specific to each lot of strip test cassettes and varies from barcode to barcode ensuring the accuracy, precision, and reliability of each test.

The lot number and test type displayed on the Vertu reader must match the label information on the strip test package. If the reader has just been turned on, confirm that “Scan” is selected on the display screen and then press the center button on the keypad. If the reader has already run a test, use the arrow keys to select “NT” and then press the center button.

When the scanner is ready to read the barcode, the reader will beep 3 times. Press the yellow trigger on the barcode scanner and aim the laser beam point at the barcode for the test you are running. When scanning the barcode, it may be necessary to adjust the distance between the scanner and the barcode while pressing the yellow trigger. When the barcode is scanned, the reader will beep once and the display screen will show the correct information.

There is a different bar code for ground peanuts and peanut paste.

PREPARATION OF PREMIXED AQUEOUS EXTRACTION SOLUTION

Water plus AQUA Solution A and AQUA Solution B can be added individually to each sample when running one sample at a time. Alternatively a premixed Aqueous solution containing all three reagents (water, AQUA Solution A and AQUA Solution B) can be made in advance and the one solution added to the sample.

To make 900 mL of AQUA Premix Solution, combine the following reagents in a glass or plastic container in the order listed below:

1. 540 mL water
2. 180 mL AQUA Solution A (one bottle VICAM p/n 100000336)
3. 180 mL AQUA Solution B (one bottle VICAM p/n 100000337)


Mix well.

Alternatively, the following AQUA Premix solutions can be purchased from VICAM:

- 900 mL (for 18 tests) p/n 100000339
- 4 L (for 75 tests) p/n 100000345
- 20 L (for 400 tests) p/n 100000347

PRODUCT FEATURES

Afla-V Strip Test Cassette

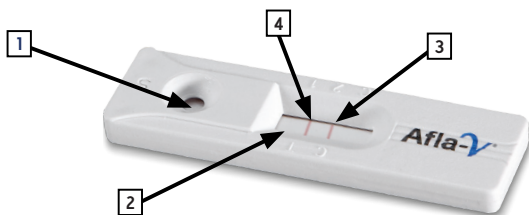

- 1 Sample Well
- 2 Results Window
- 3 Control Line
- 4 Test Line

PIPETTE TERMS

Aspirate – to draw the sample up into the pipette tip

Blow out – to empty the tip completely

PIPETTING TECHNIQUES

1. Press the operating button to the first stop.
2. Dip the tip into the solution and slowly release the operating button to "wet" the tip.
3. Press the operating button to the first stop again and slowly release the operating button. Wait 1–2 seconds and then withdraw the tip from the liquid, touching it against the edge of the reservoir to remove excess liquid.
4. Dispense the liquid into the receiving vessel by gently pressing the operating button to the first stop and then press the operating button to the second stop. This action will empty the tip. Remove the tip from the vessel, sliding it up the wall of the vessel.
5. Release the operating button to the "Ready" position.
6. Please consult the chart for visual instructions on operating the pipette. Be sure to always hold the pipette in a vertical position.

PIPETTE POSITIONS

1. Press pipette button down to first stop and place into liquid.

2. With tip in liquid, slowly aspirate (draw up) by letting button return to "Ready" position.

3. Dispense liquid into other vessel by pressing button to first stop and then to second stop. This will blow out the tip.

4. Lift the tip out of liquid, and blow out any remaining liquid by releasing and pressing the operating button again. Return button to the "Ready" position.



1. Calibrate the Vertu reader daily by scanning in the Afla-V AQUA Peanut bar code for ground peanuts or the Peanut Paste bar code for peanut paste for the lot of strips being used. You will hear a beep when the bar code is accepted. Make sure the lot number displayed on the reader is the same lot number as the strips.
2. Weigh 10 g ground sample into an extraction tube.
3. Add to the extraction tube 50 mL of AQUA Premix Solution.
(Alternatively, 30mL of purified water followed by 10 mL of AQUA Solution A, then 10 mL AQUA Solution B can be added to the sample.)
4. Cap the extraction tube and vortex at high speed for 2 minutes.
5. Filter the extract into a clean extraction tube.
6. Transfer 100 μ L of the filtered solution to the circular opening on the Afla-V strip at a rate of about 1 drop per second. Be sure to hold the pipette vertically. The pipette tip can touch the bottom of cassette.
7. Allow the strip test to develop for 5 minutes.
8. Insert the Afla-V strip test into the Vertu reader (circular opening side in first).
9. Press the center key on the reader to take a reading.
10. To print the result, use the arrow keys to move the cursor to “P” in the upper left-hand corner of the display screen. Then press the center button on the keypad.
11. To run the next sample, use the arrow keys to move the cursor to “NT” and press the center button on the keypad. If using the same lot of strip test cassettes, insert the next sample. If using a different lot, scan in the new barcode before inserting the next sample.

Assay range: 0–100 ppb

Limit of detection: 5 ppb



1. When scanning in the barcode, make sure the red + in the center of the barcode scanner is positioned in the center of the barcode. It may be necessary to adjust the height of the barcode scanner to position it correctly for scanning.
2. Press the center button on the keypad quickly and gently to read samples. Pressing the center button on the keypad long and hard will turn the instrument off.
3. Purified water can be reverse osmosis purified, distilled or deionized water.
4. When running one sample at a time, add water first, then AQUA Solution A and AQUA Solution B.
5. All strips should have a visible control line.
6. Do not run test in a location where air from an air conditioner, heater, or window will blow directly on strips.



BENEFITS

Sustainable – Safe, solvent free testing

Fast screening – Results in 5 minutes*

Simple – No special training required

Precise – Limits of detection as low as 5 ppb

Convenient – Easily performed onsite or in the lab

Durable – Long shelf life

Accurate – Real-time data which can be printed or downloaded to a computer

Wide Range – 0 to 100 ppb

**After extraction.*







A Waters Business

KEY LOCATIONS

Headquarters:

34 Maple Street
Milford, MA 01757
USA
Tel.: +1 800 338 4381
+1 508 482 4935
Fax: +1 508 482 4972

Orders:

1848 N. Deffer Drive
Nixa, MO 65714
USA
Tel.: +1 877 228 4244
+1 417 725 6588
Fax: +1 417 725 6102

www.vicam.com

Waters

THE SCIENCE OF WHAT'S POSSIBLE.®

Subject to change without notice.

©2016 Waters Corporation. Waters, The Science of What's Possible, VICAM, Afla-V and Vertu are registered trademarks of Waters Corporation. AQUA and Cultivating Success Through Science are trademarks of Waters Corporation.

The analytical methods presented in this data sheet have been researched and developed by VICAM to be used exclusively with Afla-V AQUA products. These methods have been validated in the VICAM laboratories to perform to the specifications indicated in the Afla-V AQUA procedures. The user assumes all risk in using Afla-V AQUA procedures and products. VICAM makes no warranty of any kind, expressed or implied, other than that Afla-V AQUA products conform to VICAM's printed specification and quality control standards. VICAM will, at its option, repair or replace any product, or part thereof, which proves to be defective in workmanship or material. VICAM's undertaking to repair or service such products is exclusive and is in lieu of all other warranties whether written, oral, expressed, or implied, including any implied warranty of merchantability or fitness for a particular purpose. VICAM shall have no liability for anticipated or lost profits or any loss, inconvenience or damage whether direct, indirect, incidental, consequential or otherwise, to person or property, or for strict liability or negligence arising from or in connection with the use of these assay procedures or Afla-V AQUA products.